

Many text books teach how to do various gas chromatographic configurations to allow the measurement of multiple components in several type of background. Most of the following configuration and text was adapted from the book "The analysis of gases by chromatography" by C.J. Cowper and A.J. Derosé, Pergamon Press. Another source of useful information was found in "Fundamentals of Gas Analysis by Gas chromatography" by Brian Thompson of Varian Associates Inc. As far we know both publications are out of print.

All of the following configuration could be done with diaphragm or standard rotary valves. However the best performance are achieved with our CLP type rotary valve. This is the reasons why all valve drawing show the internal purging grooves that characterize the CLP series valves. Long term column or sample loop isolation are best achieved with the MDVG or DV series diaphragm valves. Column and sample could be isolated for hours without noticeable effect. All suggestion are welcome and could be also published. We enjoy customers feedback.


## COLUMN SELECTION

### App. #1

A four-port valve is used as shown in Fig. 1. Operation of the valve allows the injected sample to go onto either column. The two detectors could be the sensor and reference cells of a thermal conductivity detector.

When used in this way the system may not offer any advantage over the use of another gas sampling valve. With a gas sampling valve in each carrier gas line, samples can be injected onto either column without the baseline disturbances which may attend the use of switching valve.

If the effluent from a single column is to be switched to different detectors during an analysis, then the configuration of Fig. 1 can be modified by fitting the column before the valve. The effluents from the two columns can then be switched between the two detectors.

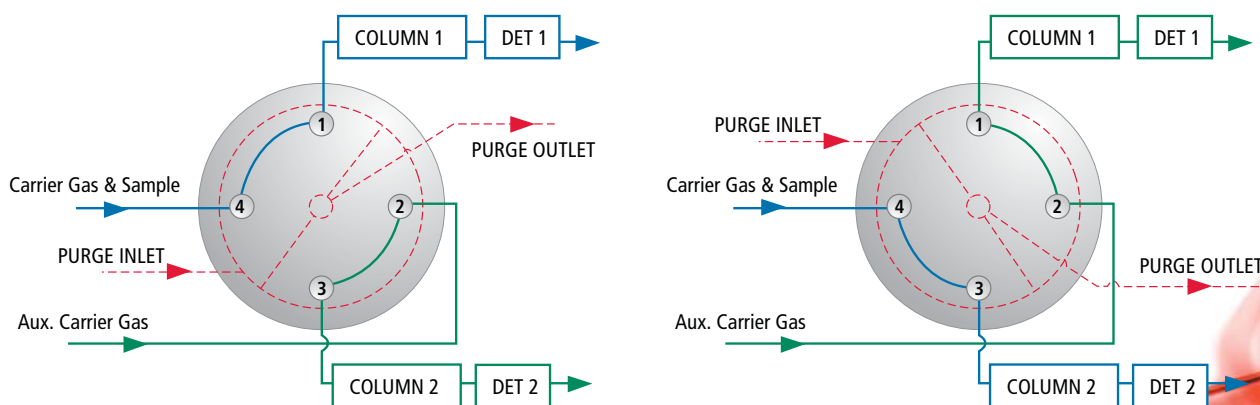

■ Fig. 1: Column selection

## BACKFLUSHING


If the sample is complex, optimising the separation of a group of components may mean that heavier components have very long retention times. Conditions may not favour their measurement and they can cause an excessive delay between analyses. The gas flow in the column can be reversed after elution of last component of interest, and the heavier components eluted backwards: this should take approximately as long as they have spent travelling forwards in the column.

The backflushed components can be vented, or fed to the detector so that a total figure for heavy components may be measured. While components are being backflushed, they are to some extent regrouped: those components which travel faster in forward flow also travel faster in backflush, and so catch up with the slower-moving ones. This is a somewhat simplified view and depends on the pressure drop and the ratio of retention times of the heavier components.

Peak broadening cannot be reversed, but the backflushed components often appear as a single, perhaps rather unsymmetrical peak. The limit of detection of this group is obviously much lower than that of the individual components which comprise it.

### Backflush column to detector

A four-port valve is used as shown in Fig. 2. If the detector is an FID, the pressure surge when flow is reversed may extinguish the flame: a restrictor should be fitted between the valve and the detector.

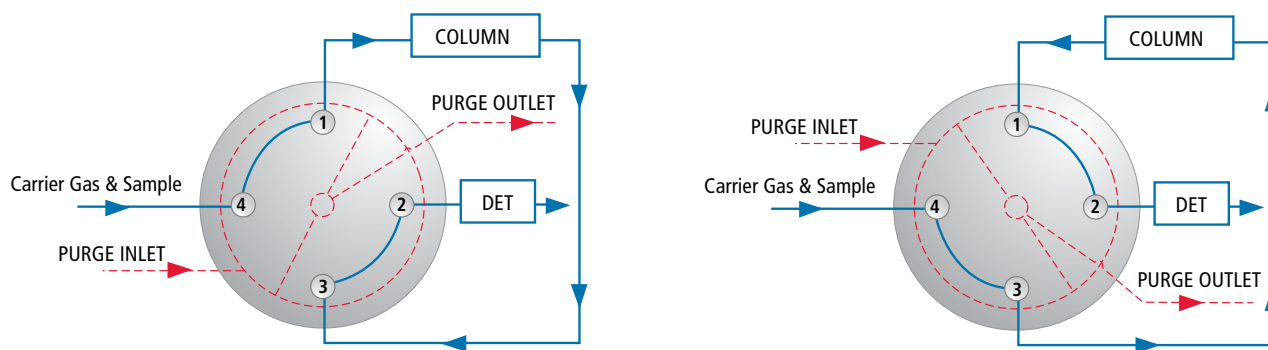


■ Fig. 2: Backflush column to detector

## Backflush pre-column

With the above configuration, the total time taken is approximately twice that of the forward flow. An alternative approach uses two columns, the first of which (pre-column) need only be long enough to separate the group of desired components from the group to be backflushed. The first group is passed to the second (analytical) column in which they are separated. Simultaneously, the second group is backflushed from the first column. A six-port valve is used as shown in Fig. 3.

Restrictors R1 and R2 are set to have similar pneumatic resistances to the analytical column and the pre-column respectively. This should ensure that the flow through the analytical column is undisturbed by the operation of the valve.

The backflushed group of components can be measured, if desired, by placing another detector on the vent line. Alternatively the vent line can be teed in to the existing detector after the analytical column. In this case the column lengths and switchings time must be chosen so as to avoid interference between the backflushed group and the normally eluting components.

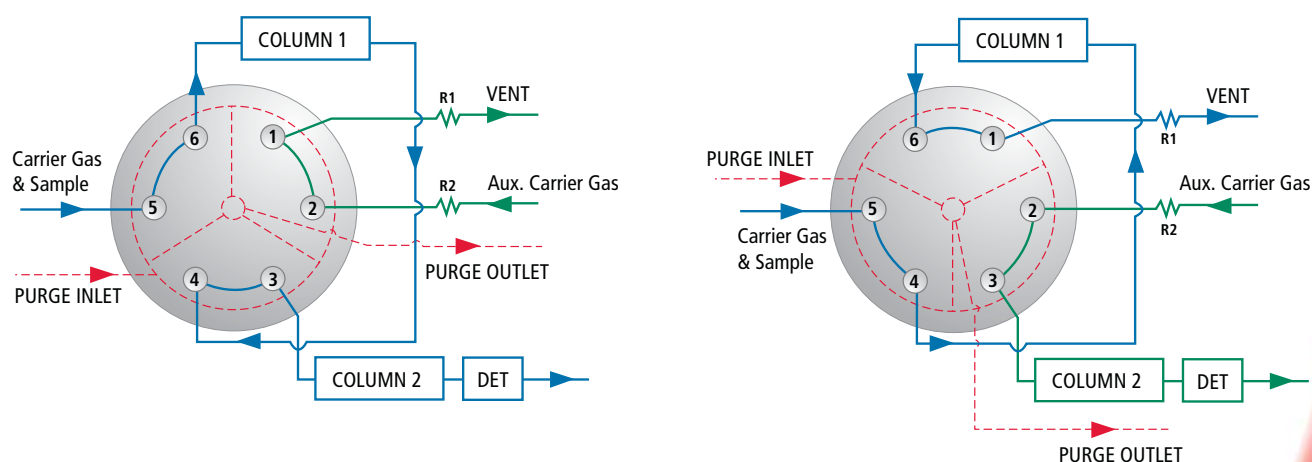


■ Fig. 3: Backflush pre-column to vent

## Backflush and sequence reverse

A six-port valve is used a pre-column and analytical column. A single carrier gas source is used, and the plumbing arranged so that switching both reverses the flow through the pre-column and positions it after the analytical column. Fig. 4 shows the configuration.

The pre-column would normally be chosen to be considerably shorter than the analytical column. This means that it has little effect on the normally eluted components, even though they pass through it twice. The backflushed group of components, however, having spent only a short time travelling in either direction will emerge as a sharp peak. It can be arranged for the backflushed group to be the first to elute: this allows much lower detection limits than with a more conventional broad backflushed peak.

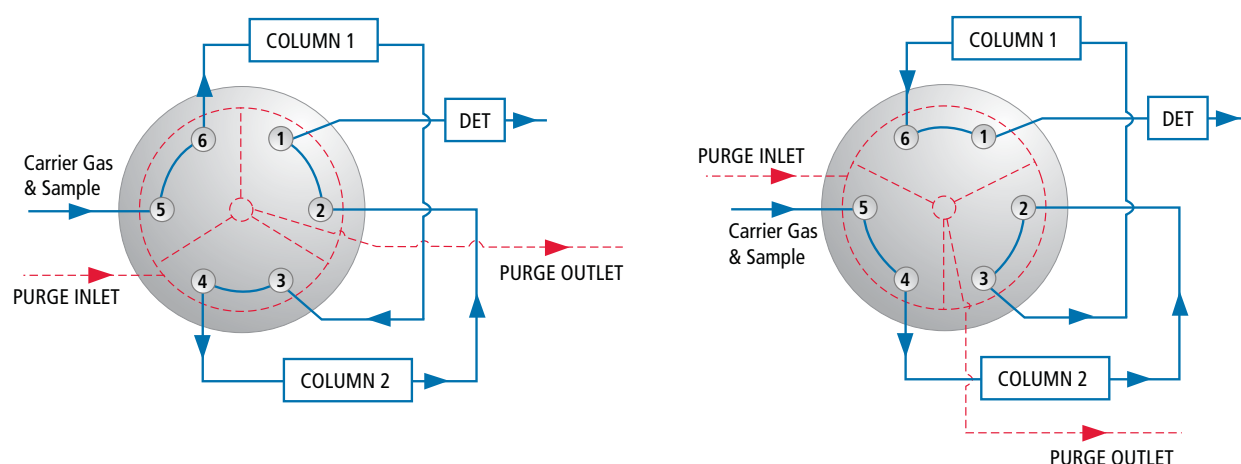


■ Fig. 4: Backflush and sequence reverse

## App. #3 COLUMN ISOLATION

Series operation of columns allows measurement of groups of components separated on columns whose characteristics are different. The drawback of the technique is that all the components eluted from the first column pass to the second. This can mean excessive retention times, which prolong the interval between analyses, or deactivation of the second column (e.g. H<sub>2</sub>O, CO<sub>2</sub> on molecular sieve).

A switching valve can be used to isolate and by-pass the second column after the appropriate components have entered it. Elution proceeds from the first column directly to the detector, then subsequently the second column is reconnected and the components which have been stored in it are passed to the detector. Alternatively, it may be possible to continue elution through the columns in series until all components from the second column have been measured, then by-pass the second column before the heavier components have reached it.

Fig. 5 shows the use of conventional six-port valve. The restrictor is adjusted to have the same resistance as the column, so that the flow is independent of the valve position. A disadvantage of the configuration shown in Fig. 5 arises from the fact that both ends of the column are connected via the valve when it is isolated. During isolation, the pressure of the carrier gas in the column will adopt a value intermediate between the inlet and outlet pressures under flow conditions. This is achieved two ways: by continuing forward flow through the column, and by reverse flow from the high-pressure to the low-pressure end of the column via the valve. There is a danger that a component which has just entered column 2 before it is isolated may be transferred after isolation to the low pressure end of the column. This is avoided by the configuration shown in Fig. 6.

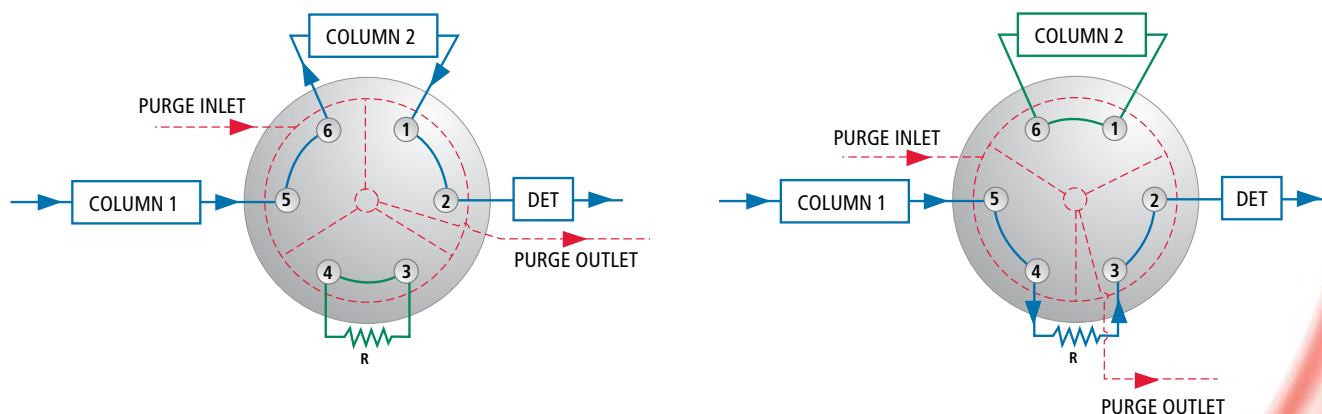


■ Fig. 5: Column isolation - method 1



## COLUMN ISOLATION (CONT'D)

Here the use of a rotor with only two slots means that equalisation of pressure in the isolated column is only achieved by forward flow of carrier.

When components are isolated in a column in this way, the chromatographic process stops, but the component bands are still subject to diffusion in the mobile phase, which causes some broadening. The degree of broadening depends upon the time and upon the partition coefficient of the component: significant band broadening only occurs in the mobile phase and hence affects lighter components to a greater extent. In general, within the time-scale of most analyses, the degree of broadening will be perceptible but not troublesome. Furthermore, this type of operation will almost invariably be controlled by an automatic timer, and so the peak shape and resolution will be consistent.

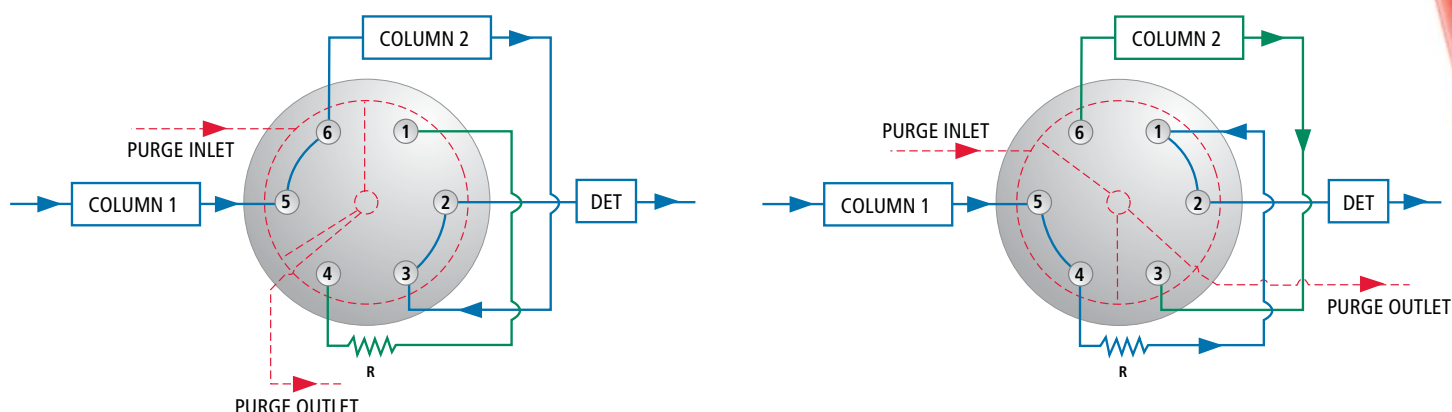


■ Fig. 6: Column isolation - method 2

## FOREFLUSHING

Column isolation allows maximum flexibility when multiple columns are being used, as components can be stored until the most appropriate time in the analytical cycle. There is the danger, however, that if the isolating valve is not completely leak-tight, stored components may leak out or small amount of air may leak in. The CLP series eliminate these problems. Foreflushing has been described by Willis (1978) as a way of avoiding this. Fig. 7 shows the use of a six-port valve for this purpose.

After sample injection, and when the lighter components have been passed on to column 2, the valve is rotated, and the heavier components eluted directly from the column 1 to the detector. The lighter components, having been separated on column 2, now pass through column 1 for a second time and are then detected. Provided that they have been sufficiently well separated on column 2, the small extra amount of peak broadening caused by this second passage through column 1 is not significant. Column lengths must be selected so that there is no overlap of components.

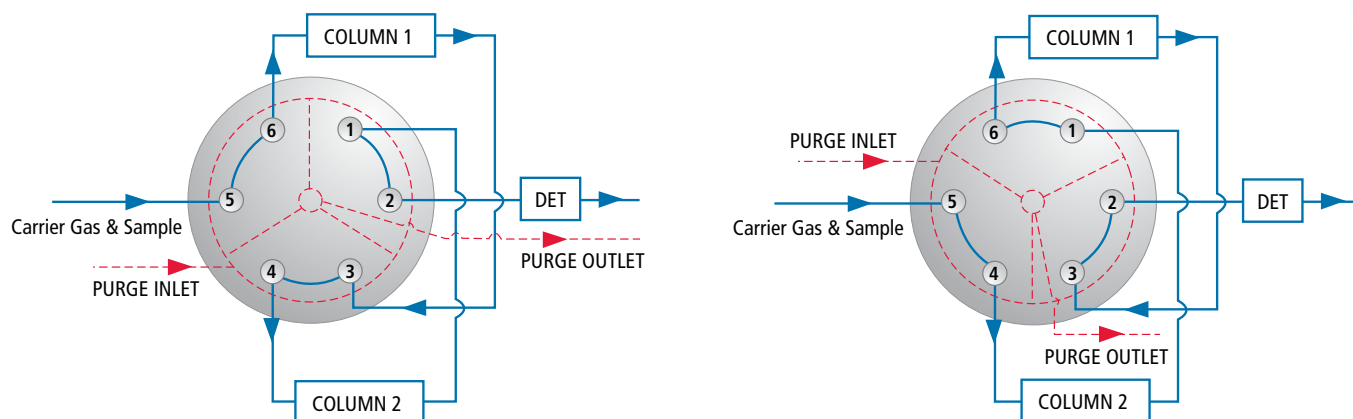

■ Fig. 7: Foreflushing


## MULTIPLE FUNCTIONS

When an analysis requires more than one of the above operations, or a combination of any of these with sample injection, a series of valves can be used. Alternatively where two functions can occur at the same time, they can be dealt with by a single multi-port valve. Some typical examples follow.

### Gas sampling and backflush to detector

An eight-port valve with 45° rotation is used as shown in Fig. 8. The left hand position shows the sample being loaded into the loop. The right hand position shows the sample being injected and flowing through the column in what would be considered to be the forward direction. Components separated in this way continue on the further columns or to the detector. When the valve is returned to the left hand position the column is backflushed and the sample loop returned to the load position.

The timing of the backflush operation is dictated by the requirements of the analysis. The time at which the sample loop is returned to the load position is not critical, and so the fact that these operations necessarily occur simultaneously is not disadvantage.

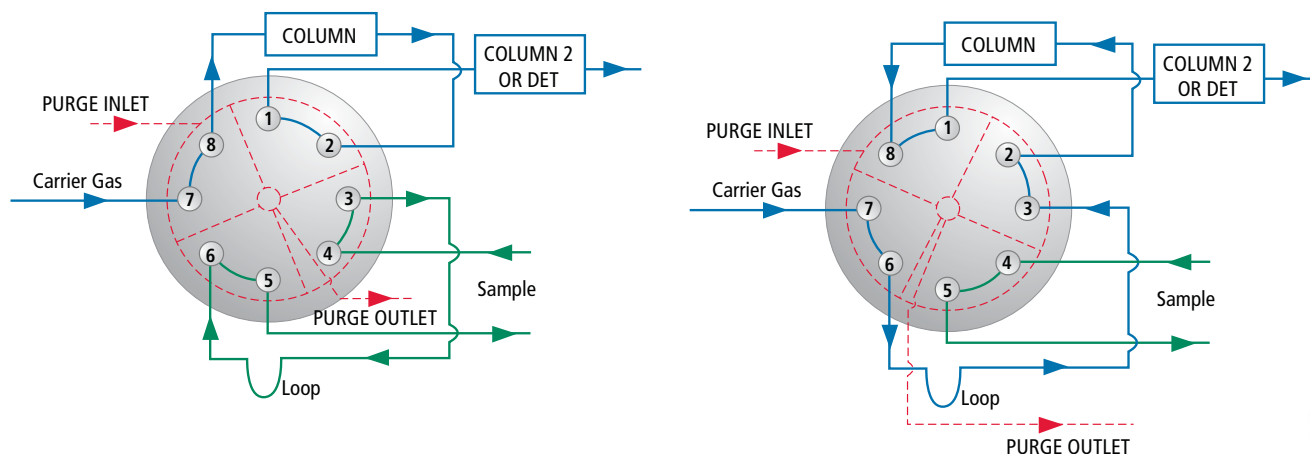


■ Fig. 8: Gas sampling and backflush to detector



## Gas sampling and backflush to vent


A ten-port valve is used with an auxiliary supply of carrier gas as shown in Fig. 9. Restrictor 1 is adjusted to have the same resistance as Column 1, and restrictor 2 to the same resistance as any downstream columns.

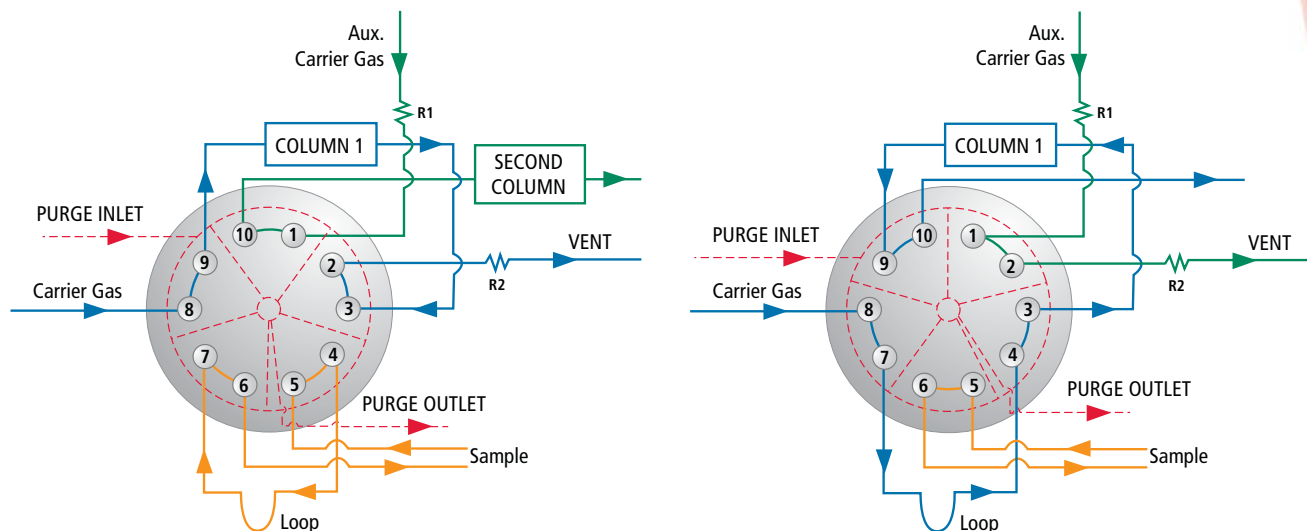

■ Fig. 9: Gas sampling and backflush to vent

## Gas sampling and sequence reverse


A ten-port valve combines these functions as show in Fig. 10.

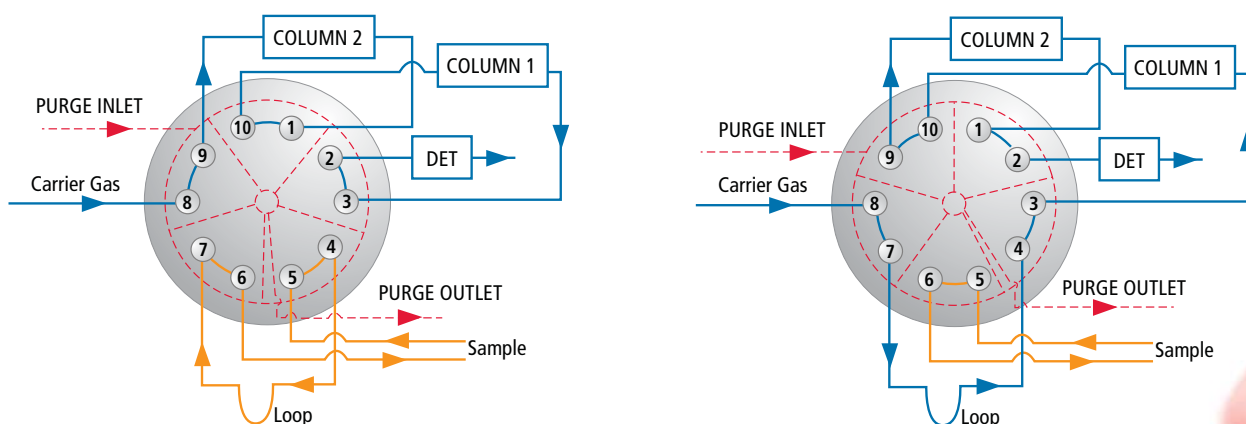

■ Fig. 10: Gas sampling and sequence reverse

### Column selection and backflush

Either of two columns is selected: the other is simultaneously backflushed. An eight-port valve with 90° rotation is used as shown in Fig.11.

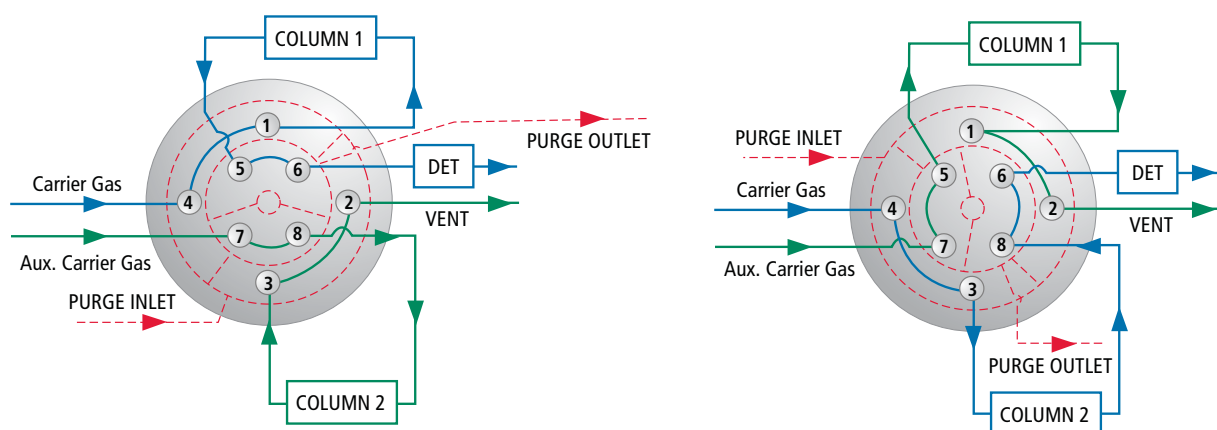


■ Fig. 11: Column selection and backflush

## GROUP ISOLATION / ENHANCE SELECTIVITY

This configuration use two different type of column and two different size of sampling loop. Normally SL1 is smaller than SL2. This allow to transfer a poorly resolved peaks from column 1 to SL2, and then re-inject SL2 into column2 that have appropriate separation for the peaks of interest.


This work better than standard heartcut since SL2 is quickly injected into the column 2 reducing peak broadening and improving separation. It may also reduce the load on column 2 by first pressurizing SL2 before to inject it.

This is simply a matter of timing. Optionally an auxiliary detector could be added on V2 vent line for parallel measurement of impurities resolved by column 1. The addition of Valve 4 allow fast tuning of the system, this could be useful for LAB GC where configuration are frequently change.

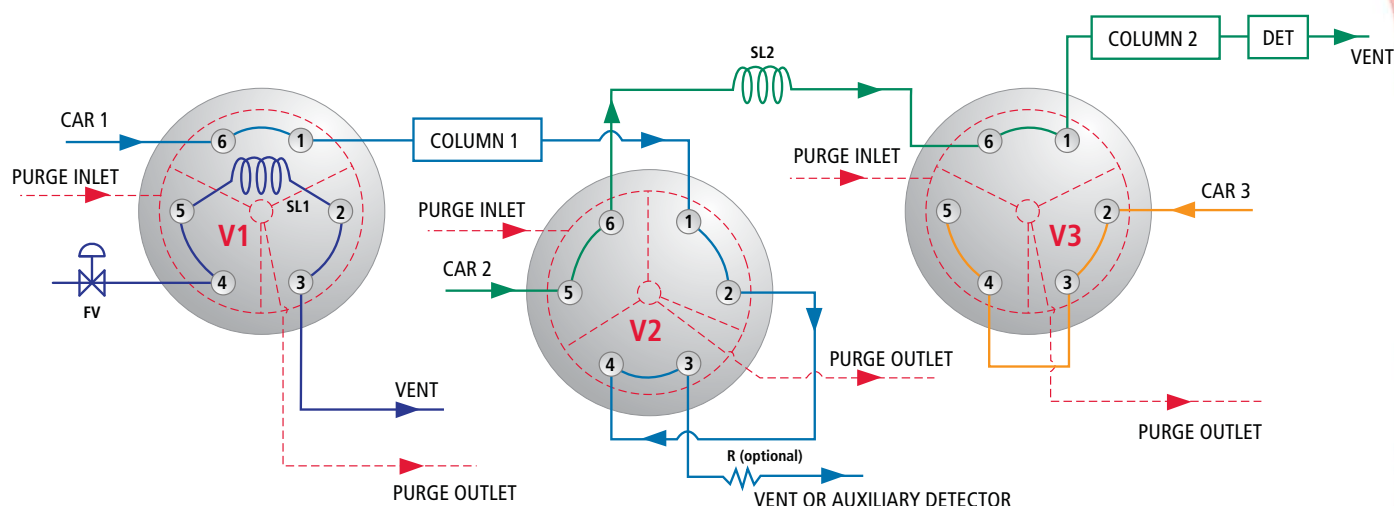


■ Fig. 12: - Filling SL1  
- Injecting SL2 on column 2  
- V3 port #5 plugged

## GROUP ISOLATION / ENHANCE SELECTIVITY (CONT'D)

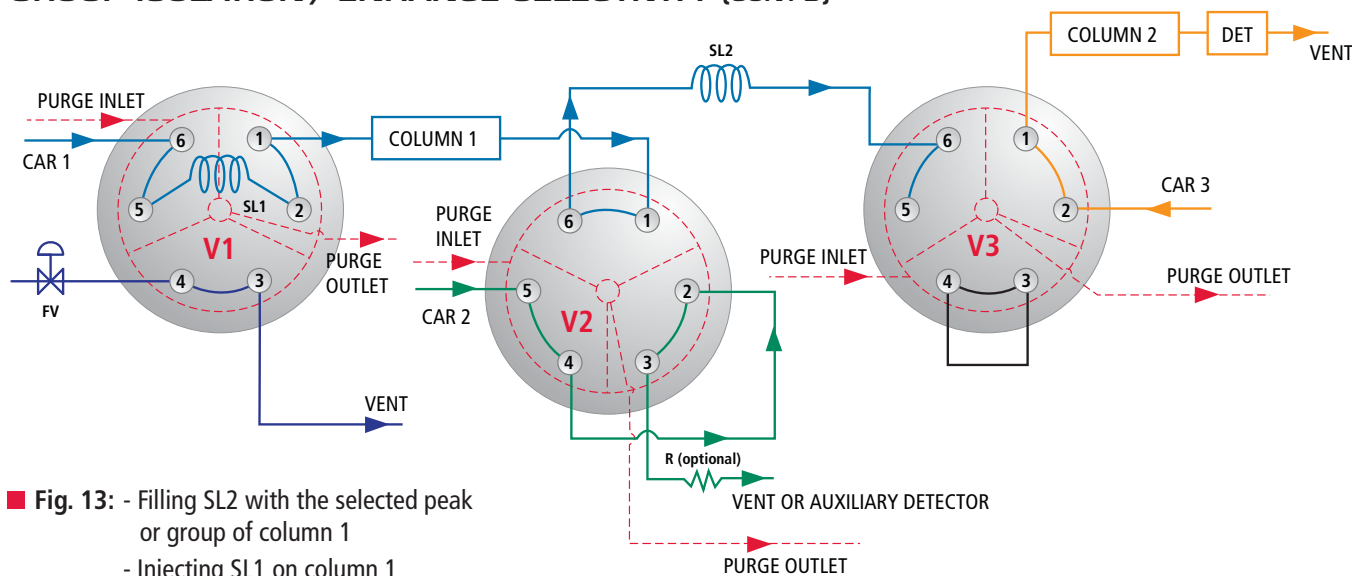

■ **Fig. 13:** - Filling SL2 with the selected peak or group of column 1  
- Injecting SL1 on column 1

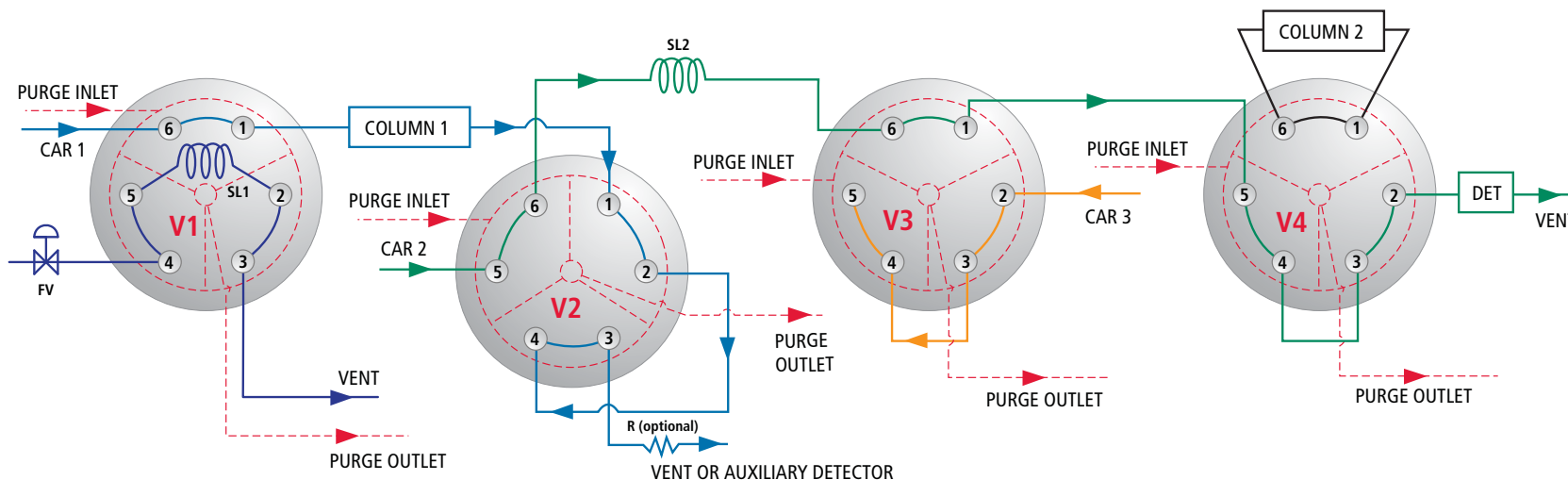

■ Fig. 14: Configuration variation

Here V4 is added to allow column 2 by-pass. This is a useful tool to allow proper timing FV3, in order to synchronize V3 operation.

## DISCLAIMER

Although Analytical Flow Products ("AFP") endeavours to ensure the accuracy of the information provided within the present Application Notes (the "Notes"), said Notes are provided strictly on an "as is" basis and no representation, warranty or guarantee, express or implied, is made or given as to (i) the accuracy or appropriateness of said Notes for any given purpose or circumstance (ii) whether or not the information or material displayed within the Notes may or may not infringe the rights of any third parties or (iii) whether any trade name or trade mark associated to third party products and mentioned within the Notes belong to the owner of the mark or name.

It is the reader's responsibility to analyse, evaluate and validate the information provided in the Notes in a responsible and professional manner before deciding whether or not to use said information in any way and thereby assume all risks associated with such decision.

AFP assumes no liability or responsibility whatsoever with respect to (i) any errors or omissions which may be present within the Notes, (ii) any infringement of third party rights, including patents or other intellectual property rights arising from the use of the information contained within the Notes, or (iii) any loss or damage (be it direct, indirect, consequential or otherwise) whatsoever which may in any way arise out of or in connection with the use or application of any information provided within the Notes, including, without limitation, any errors, falsehoods, or inaccuracies which may be contained therein.

Unless otherwise specified, AFP is the owner of copyright in the Notes and same cannot be used or reproduced in any way without obtaining AFP's prior written consent. AFP reserves the right to modify, update or deny access to the Notes at any time.

