LYOMAC FREEZE DRYER

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Freeze dryer series

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Level analysis experimentation congealer series

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Zhejiang Lyomac pharmaceutical equipment Co.,LTD. was founded in 2009 and specializes in the R&D, production and sales of vacuum freeze-drying products. With the vacuum freeze dryer as the core product, forming different scales and different applications of three series of products as well as targeted solution for scientific research experiment, medical, and food industry. We are well-known on providing the vacuum freeze-drying equipment and targeted solution in China. We are one of the first domestic enterprises engaged in vacuum freeze-drying technology and safety production standardization of Beijing industrial enterprises. Our head office is in Shanghai. We have the branches and special maintenance stations in Shanghai, Guangdong, Wuhan, Xian, Chengdu, and Hainan. Our products are exported to the United States, Russia, South Korea, Japan, Singapore, Malaysia, Canada, Brazil, Argentina and other countries.

The company has developed three major series of products with independent intellectual property rights and high-tech content. In 2012, the company researched one kind of patented shelf for vacuum freeze dryers, which is widely used in the current mainstream pilot freeze dryers and production freeze dryers. It makes the rapid development of domestic vacuum freeze-drying technology. In 2015, we researched and developed a kind of energy-saving vacuum freeze-dryer system and a vacuum freeze dryer with this system. At present, this energy-saving system is mainly used in big scale freeze-drying equipment, effectively reducing energy consumption of equipment by about 30%.

With Kind, open, rigorous corporate culture and scientific management, Songyuan Huaxing has excellent professional managers and management team. Senior refrigeration experts, senior vacuum experts, senior electronic control experts and senior mechanical design experts form technical teams. They cooperate close in the product development, promotion and maintenance. They have a profound understanding the meaning of "customer-centric". "service" is carried out throughout operation and management, to win wide trust and support from customers. With an open and harmonious attitude, Songyuan Huaxing has a deep technical cooperation with many well-known freeze-drying technology manufacturers and universities at home and abroad.

The process of growing up is the process of serving customers and making success with the customers together. We try our best to do everything well and meet every challenge with confidence.

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Vacuum freeze dryer is widely used in medicine, pharmacy, biology research, chemical industry and food production, etc. After freeze drying process, a long term preservation for material is much easier. They can be restored to original state and maintain their chemical and biological characteristics after being watered.

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LGJ-10/12/18 freeze dryers, small amount of samples can pre-freeze in condenser. They are suitable for lab use or small-scale production, and they can meet requirements for most laboratories.

Types : A : standard type ,B: top press type, C: manifold type , D: manifold with top press type, E :T-type

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Main features:

- All-in-one structure, small size, no external flange, easy to use, no leakage.
- All materials in contact with the product use inert materials to meet the requirements of the GLP.
- The cold trap and console are made of stainless steel, which is anti-corrosion and easy to clean.
- All stainless steel inflated/water release valve is designed and produced by our company, which is safety, anti-corrosion, no leakage.
- With sample pre-frozen function, large cold trap opening, no inner coil, and low temperature refrigerator is needless.
- Patented gas diversion technology, and strong ice holding ability.
- World - known brand compressors, high efficiency, long life, low noise.
- Well-known brand vacuum pump with a high pumping speed to achieve a higher ultimate vacuum.
- Vacuum pump protection function can set the cold trap temperature to protect the vacuum pump life.
- Professionally designed FD-LAB freeze-drying machine control system + SH-HPSC-I modular controller with high reliability and stability.
- Intelligent data recording system can real-time record and display the cold trap temperature curve, sample temperature curve, vacuum curve.

预冻功能
Pre-freezing
function

思科普压缩机
Secop compressor

混合环保制冷剂
Green environmental
protection compound
refrigerant

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Model Selection Guide:

- A Standard:
It is suitable for freeze-drying bulk (liquid, pasty, solid) conventional materials.
- B Top-press:
It is not only suitable for freeze-drying of bulk (liquid, pasty, solid) conventional substances, but also suitable for drying of Vials materials. When preparing for freeze-drying, the materials are separated into Vials according to the need. After floating the cap, the freeze-drying process start. After drying, the capping device is operated to tighten the cap, which can avoid secondary pollution, re-adsorb water, and is easy to store for a long time.
- C Multi manifold general type:
It is suitable for the freeze-drying of bulk (liquid, pasty, solid) conventional materials, and can be fitted with flasks outside the drying chamber to dry the materials rotated inside the bottle wall. At this time, the flask is connected to the manifold outside the drying chamber as a container. The material in the flask is heated at room temperature. Through the multi-manifold switch device, the flask can be removed or loaded at any time as needed without stopping the machine.
- D Multi manifold cover type:
On the basis of the characteristics of the common type, it combines the characteristics of the capping type and the common type of multi-weir pipe. It is suitable for freeze-drying of bulk (liquid, pasty, solid) conventional material ; suitable for drying of vials materials. When preparing for freeze-drying, the materials are separated into vials according to the need. After floating the cap, the freeze-drying process start . After drying, the capping device is operated to tighten the cap, which can avoid secondary pollution, re-adsorb water, and is easy to be stored for a long time. The flask is used to dry the material whirled inside the flask. At this time, the flask is connected to the manifold outside the drying chamber as a container. The material in the flask is heated at room temperature. Through the multi-manifold switch device, the flask can be removed or loaded at any time as needed without stopping the machine.
- E T shelf :
It is suitable for the preservation of microorganisms such as bacteria and fungi in ampoule tubes. The ampoule storage method is often used in laboratories. The samples were packaged in small ampoules after freeze-drying. It has the advantages of Long-term preservation, small variation, easy storage and Widely used.

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型号 Model		LGJ-10				LGJ-12/12S		LGJ-18/18S	
规格 Specification	单位 Unit	普通型 Standard	压盖型 Top-press	普通型 Standard	压盖型 Top-press	普通型 Standard		压盖型 Top-press	
冻干面积 Freeze-drying area	m ²	0.12	0.08	0.12	0.08	0.18	0.27	0.09	
冷阱温度 Cold trap temperature	℃	< -56 (空载) , 可选 < -80 (空载) < -56 (no-loading), option < -80 (no-loading)							
真空度 Vacuum degree	Pa	≤ -56 (空载) ≤ 5 (no-loading)							
捕水能力 Water-capture capability	kg/24h	3-4			3-4		6		
物料盘尺寸 Tray dimension	mm	φ 200	φ 180	φ 200	φ 180	φ 240	φ 240	φ 200	
物料盘数量 Trays number	个	4	3	4	3	4	6	3	
物料盘间距 Tray spacing	mm	70			70		70		
整机功率 Power	w	970			970		1400		
主机重量 Weight	kg	41			62		105		
主机外形尺寸 (-56℃) Dimension (-56℃)	mm	615 × 450 × 370			580 × 500 × 720		630 × 580 × 970		
主机外形尺寸 (-80℃) Dimension (-80℃)	mm	850 × 680 × 405			770 × 550 × 720		810 × 580 × 950		
干燥室尺寸 Drying chamber size	mm	Φ260 × 430	Φ260 × 490	Φ260 × 430	Φ260 × 490	Φ300 × 445		Φ300 × 540	
冷阱直径 Diameter of cold trap	mm	Φ215			Φ215		Φ270		
冷阱深度 Cold trap depth	mm	140			240		400		
可装物料 (料厚 10mm) Bulk capacity (thickness 10mm)	L	1.2	0.8	1.2	0.8	1.8	2.7	0.9	
可装入西林瓶 Vial loading number	支	—	Φ12mm: 492	—	Φ12mm: 492	—	—	Φ12mm: 615	
	支	—	Φ16mm: 297	—	Φ16mm: 297	—	—	Φ16mm: 345	
	支	—	Φ22mm: 147	—	Φ22mm: 147	—	—	Φ22mm: 183	
除霜模式 Defrosting	—	自然化霜 (可选电除霜) Natural defrosting (option electric defrosting)							
电源 Power	—	单相 3 线, 220 伏 1Ph + N + PE, 220V							

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LGJ-12S/18S freeze-dryer is a small vertical freeze-drying equipment, which has the functions of baffle heating and programmable. They are suitable for lab use or small-scale production, and they can meet requirements for most laboratories.

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Main features:

- Integral structure design, small size, no external flange, easy to use, no leakage.
- All materials in contact with the product are inert materials to meet the requirements of GLP.
- Cold trap and operating table are made of stainless steel, which is corrosion-resistant and easy to clean.
- All stainless steel inflatable (drainage) valves designed and manufactured by our company are safe, anticorrosive and leak-free.
- The cold trap has large opening, no inner coil, with sample pre-freezing function, and no cryogenic refrigerator is needed.
- Specialized gas diversion technology, uniform cold trap ice capture, ice capture ability.
- International well-known brand compressor, high efficiency and energy saving, long service life, low noise.
- Well-known brand vacuum pump, high pumping speed, achieve higher vacuum limit.
- Vacuum pump protection function, can set the starting cold trap temperature of vacuum pump, protect the service life of vacuum pump.
- 7-inch real-color industry embedded touch screen + SH-HPSC-II modular controller, high control accuracy, stable and reliable performance.
- Intelligent data recording system can real-time record and display the cold trap temperature curve, sample temperature curve, vacuum curve.
- The shelf electric heating safety.The control systemvoltage (36V) temperature can control the heating rate.
- Flexible manual + automatic control mode, manual groping process, automatic batch production.
- Real-time alarm display and historical alarm inquiry functions facilitate fault diagnosis and equipment maintenance.
- User level and password can be set, and operation management can be decentralized.
- Powerful sensor calibration function ensures long-term use of measurement accuracy.

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LGJ-10FD/30FD/50FD electric-heating freeze dryer

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LGJ-10FD/LGJ-30FD/LGJ-50FD series of freeze dryers with electric heating function (patent machine). Products can be freeze dried in situ. It changes the complex operation and prevent sample pollution, and make drying automation. These models has shelf heating and programmable functions. They can records freeze drying curves, data can be output by U-disk. User can clearly view the freeze drying process.

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Main features:

- The company's patented product square warehouse in situ vacuum freeze dryer, pre-freezing, drying in situ completed, easy to operate, good drying effect.
- The drying chamber door is made of ACRYLIC material, which is colorless, transparent and can observe the whole process of freeze-drying.
- The inflatable (drainage) valve adopts safety diaphragm valve, which can connect inert gas source, fill inert gas after drying, and prolong the shelf life of materials.
- Specialized gas diversion technology, uniform cold trap ice capture, ice capture ability.
- International well-known brand compressor, high efficiency and energy saving, long service life, low noise.
- The patented technology of partition and the fuzzy control technology of partition temperature have the advantages of small temperature difference of shelf, high precision of temperature control and even drying effect.
- Powerful sensor calibration function ensures long-term use of measurement accuracy.
- 7-inch real-color industry embedded touch screen + SH-HPSC-II modular controller, high control accuracy, stable and reliable performance.
- The specially designed FD-EMB control system can preserve hundreds of formulations, each of which contains 50 temperature control sections, so as to improve the process optimization rate.
- Intelligent data recording system can record and display the temperature curve of cold trap, sample temperature curve and vacuum curve in real time. The derived data can be browsed and printed by computer and operated in a variety of ways to facilitate process optimization and drying effect verification.
- Flexible manual + automatic control mode, manual for exploring process, automatic for batch production.
- Real-time alarm display and historical alarm inquiry functions facilitate fault diagnosis and equipment maintenance.
- User level and password can be set, and operation management can be decentralized.

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Optional accessories and functions :

- Cold trap temperature -80℃
- Eutectic point tester
- Vacuum pump anti-return device
- Cold trap electric heating defrosting
- Oil mist filtration system
- Vacuum regulating system
- LYO-MEGA Software System
- Mobile App Monitoring System
- Backfill filter
- Eggplant bottle device
- International brand vacuum pump
- Non-standard customization

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LGJ-20F/30F/50F/100F/200F silicon oil heating freeze dryer

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LGJ-20F/30F/50F/100F/200F (silicon oil heating) freeze dryer as our patent machine, freezing and drying process made in situ. It changes the complex operation and prevent sample pollution,and make drying automation. These models has shelf heating and programmable functions.They can records freeze drying curves, data can be output by U-disk. User can clearly view the freeze drying process.

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
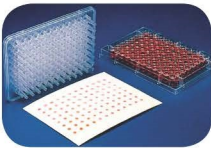
Main features :

- The company's patented product square warehouse in situ vacuum freeze dryer, pre-freezing, drying in situ completed, easy to operate, good drying effect.
- The drying chamber door is made of ACRYLIC material, which is colorless, transparent and can observe the whole process of freeze-drying.
- The inflatable (drainage) valve adopts safety diaphragm valve, which can connect inert gas source, fill inert gas after drying, and prolong the shelf life of materials.
- Specialized gas diversion technology, uniform cold trap ice capture, ice capture ability.
- International well-known brand compressor, high efficiency and energy saving, long service life, low noise.
- The patented technology of partition and the fuzzy control technology of partition temperature have the advantages of small temperature difference of shelf, high precision of temperature control and even drying effect.
- Powerful sensor calibration function ensures long-term use of measurement accuracy.
- 7-inch real-color industry embedded touch screen + SH-HPSC-II modular controller, high control accuracy, stable and reliable performance.
- The specially designed FD-PILOT control system can preserve hundreds of formulations, each of which contains 50 temperature control sections, so as to improve the process optimization rate.
- Intelligent data recording system can record and display the temperature curve of cold trap, sample temperature curve and vacuum curve in real time. The derived data can be browsed and printed by computer and operated in a variety of ways to facilitate process optimization and drying effect verification.
- Flexible manual + automatic control mode, manual for exploring process, automatic for batch production.
- Real-time alarm display and historical alarm inquiry functions facilitate fault diagnosis and equipment maintenance.
- User level and password can be set, and operation management can be decentralized.

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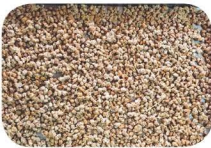Optional accessories and functions :

- Judgment function of freeze-drying end point
- Eutectic point tester
- Vacuum pump anti-return device
- Cold trap electric heating defrosting
- Oil mist filtration system
- Vacuum regulating system
- LYO-MEGA Software System
- Mobile App Monitoring System
- Backfill filter
- Data recorder
- Automatic backfill and drainage control system
- Capacitive vacuum gauge
- Eggplant bottle device
- International brand vacuum pump
- Non-standard customization

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Freeze drying laboratory：

The laboratory has a differential scanning calorimeter, scanning electron microscope, freeze-drying microscope, the above instruments for the freeze-drying process research and development provides a strong hardware support. At the same time, Songyuan Huaxing has a strong R&D team, which can provide complete technical service solutions in the following areas:

- Development of freeze-drying process
- Freeze drying process optimization
- Amplification of freeze drying process
- Freeze drying machine GMP verification

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Verification document：

Design Verification (DQ), Installation Verification (IQ), Operation Verification (OQ), Performance Verification (PQ), Factory Acceptance Test (FAT), Site Acceptance Test (SAT), Operation Manual (OM), Maintenance Manual (MM), Installation Manual (IM), Process Flow Chart (P&ID), etc..

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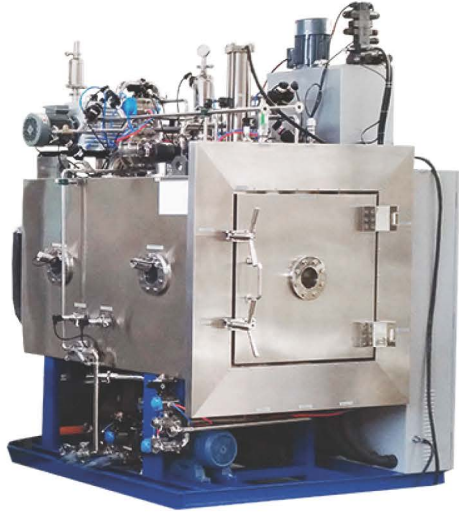
GZL/GZLY series vacuum freeze dryers are widely used in blood products, vaccines, biological products, chemicals and other pharmaceutical fields, as well as in biological research, chemical industry and food fields. Especially for Western medicine, freeze-drying machines are used in many large pharmaceutical factories. For injection, freeze-drying technology can improve the quality and storage life of medicines.

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▲ GZLY 医药型 (CIP+SIP)

Main features:

- Separators are made with proprietary patented technology with high pressure resistance, high smoothness and good temperature uniformity.
- The optimum control technology of freeze-drying curve can control the cooling rate in the pre-freezing stage and the heating rate and the vacuum value in the current stage in the sublimation and analytical drying stage.
- Professional gas diversion design and control technology, strong water catching capacity, high drying efficiency.
- The unique double-way sealing technology of mobile + fixed double-way box door sealing device sterilizes fixed silicone rubber at the same time of steam sterilization.
- Vacuum adjustment is carried out during drying operation to avoid foaming and blowing of special substances and improve drying efficiency.
- Industrial embedded touch screen + special SH-HPSC-IV modular controller, the system is stable and reliable, and the control accuracy is high.
- The specially designed LYO-CONTROL control system can preserve many sets of process formulations, and adjust the process in real time during the drying process to improve the process optimization rate.
- Flexible manual + automatic control mode, manual for exploring process, automatic for batch production.
- Customer level and password can be set, and operation management can be decentralized to meet GMP requirements.

型号 Model	单位 Unit	GZLY-1 医药型 Pharmaceutical	GZLY-2 医药型 Pharmaceutical	GZL-1 通用型 General	GZL-2 通用型 General
冻干面积 Freeze-drying area	m ²	1.15	2.16	1.15	2.16
进料量 (料厚 10mm) loading capacity (10mm thickness)	L	10	20	10	20
隔板数量 Number	层	4+1	4+1	4+1	4+1
隔板尺寸 Shelf size	mm	600 × 480	600 × 900	600 × 480	600 × 900
隔板间距 Shelf spacing	mm	80	80	80	80
控温范围 Temp.control range	℃	-50~+70	-50~+70	-50~+80	-50~+80
水汽凝结量 Ice condensing capacity	kg/B	≥ 20	≥ 40	≥ 20	≥ 40
捕水库温度 Cold trap temp	℃	≤ -75	≤ -75	≤ -75	≤ -75
降温方式 Method	-	水冷	水冷	水冷	水冷
除霜方式 Defrosting method	-	水浸泡	水浸泡	水浸泡	水浸泡
极限真空度 Final vacuum degree	pa	≤ 1	≤ 1	≤ 1	≤ 1
装机功率 installation power	kw	18	21	15	18
外形尺寸 Dimension (L*W*H)	mm	2500*1100*2000	2500*1300*2000	2500*1100*2000	2500*1300*2000
Φ22 西林瓶装量 Vial loading number	支	2460	4608	-	-
Φ16 西林瓶装量 Vial loading number	支	4788	8960	-	-
整机重量 Weight	kg	1500	2600	1200	2000
纯蒸汽耗量 SIP Pure steam consumption	kg/h	60	80	-	-
注射水量 / 纯水 CIP Water consumption	l/min	60	75	-	-

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Series of Level analysis experimentation congealer

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SL-II, SL-III level analysis experimentation congealer is special purpose low-temp congealer which used to make some experiments of level analysis, also used to other experiments in the lou-temp environment, or used to refrigerate something.

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Main Characteristics:

- Imported compressor, the refrigeration quantity is big.
- Refrigeration unit designed at machine bottom: meeting compressor's working principle, improving cooling effect, longer working life, high temperature control accuracy.
- Adopting industrial-grade true-color LCD control system, it is easy to operate and powerful.
- Professional development of SL-LAB control system, real-time display of system parameters, high control accuracy.
- The top is 1.7 meters, convenient for operation.
- All stainless steel wall, bright and anti-corrosive.
- 2 level analysis pole fixed staff, 2 levels of bearing splint.
- Own lighting lamps, antisepsis light, inner power socket.
- Own temperature exceeding, temperature shortage and alarm function.
- There is caster at bottom, easy to move.

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Technical indices:

Temperature scope	1℃ -10℃ (The difference between the upper and lower limits is not less than 4℃ .)
Temperature test accuracy	± 0.5℃
Temperature fluctuation scope	± 2.5℃
Temperature decrease time	≤ 120min.
The specification of tray	605 × 700mm (2 layers, the position of top and bottom is adjustable)
Inner cubage	800L
Power	500W
Inner dimension	660x760x1700mm
Outer dimension	760x850x1950mm

Single-door Chromatography experiment refrigerator : inner wall stainless steel, inner and shell stainless steel

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Main Characteristics:

- Imported compressor, the refrigeration quantity is big.
- Refrigeration unit designed at machine bottom: meeting compressor's working principle, improving cooling effect, longer working life, high temperature control accuracy.
- Adopting industrial-grade true-color LCD control system, it is easy to operate and powerful.
- Professional development of SL-LAB control system, real-time display of system parameters, high control accuracy.
- The top is 1.5 meters, convenient for operation.
- All stainless steel wall, bright and anti-corrosive.
- 2 level analysis pole fixed staff, 2 levels of bearing splint.
- Own lighting lamps, antisepsis light, inner power socket.
- Own temperature exceeding, temperature shortage and alarm function.
- There is caster at bottom, easy to move.

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Technical indices:

Temperature scope	1℃ -10℃ (The difference between the upper and lower limits is not less than 4℃ .)
Temperature test accuracy	± 0.5℃
Temperature fluctuation scope	± 2.5℃
Temperature decrease time	≤ 120min.
The specification of tray	505 × 615mm (2 layers, the position of top and bottom is adjustable)
Inner cubage	1200L
Power	1100W
Inner dimension	660x1200x1500mm
Outer dimension	760x1315x2000mm

Double-door Chromatography experiment refrigerator : inner wall stainless steel, inner and shell stainless steel